

# Corning® Stripettor® Pro Pipet Controller

CORNING

## Instruction Manual

Catalog Numbers:  
4999  
4900



UK  
CA CE



## Table of Contents

1.0 Introduction .....	1
2.0 Safety Information .....	2
3.0 Limitations of Use .....	3
4.0 Switching On .....	3
5.0 Aspirating and Dispensing Liquids.....	4
6.0 Troubleshooting.....	6
7.0 Replacing the Filter .....	7
8.0 Charging the Battery .....	8
9.0 Replacing the Battery .....	9
10.0 Maintenance.....	10
11.0 Components.....	10
12.0 Spare Parts.....	11
13.0 Limited Warranty.....	11
14.0 Equipment Disposal.....	12


## 1.0 Introduction

The pipet controller is a device intended for pipetting liquids with the use of measuring pipets. It can work with all types of glass or plastic pipets in the volume range from 0.5 mL to 100 mL.


Two device speed modes (LOW and HIGH) which can be chosen from 3 predefined values for each of them allow user to customize the device to the application. Two dispense modes permit selection of dispensing intensity depending on the user's needs. The selected setting of the pipet controller mode is shown on the display.

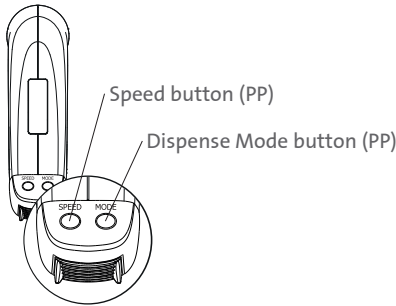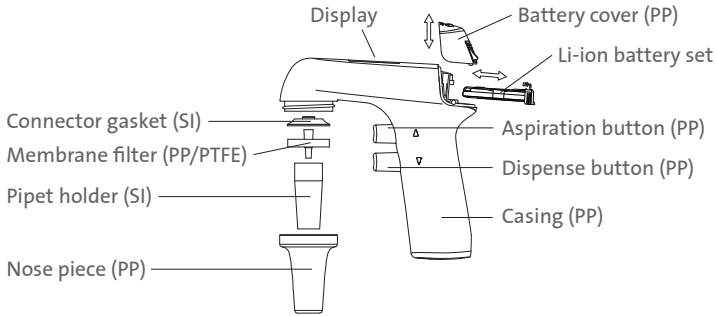


PP: Polypropylene  
PTFE: Polytetrafluoroethylene  
SI: Silicone

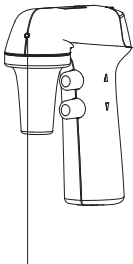

Direct outlet of liquid vapors

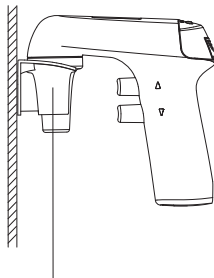

Wall hanger (PP)

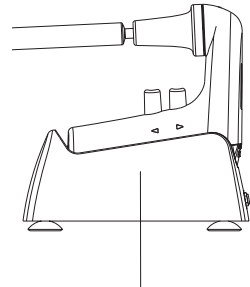

Charging stand (PP)  
(included with Cat. No. 4999)

## 2.0 Safety Information



**WARNING:** Risk of injury

**CAUTION** Risk of damage to the device or errors in pipetting of liquids.

Before starting the work with the pipet controller every user should read these operating instructions carefully.

### **CAUTION:**

- Using the device inconsistently with the operating instructions may result in damaging the device.
- The device should be serviced only at an authorized service center, otherwise the manufacturer will be relieved from any liability under the warranty.
- Only original spare parts and accessories, recommended by the manufacturer, shall be used.
- Only the original charger, supplied by the manufacturer, shall be used for charging the battery.
- In case of incorrect functioning of the pipet controller, work shall be stopped. The device shall be cleaned according to Section 9 and sent for repair to an authorized service center.
- In the case of mechanical damage to the casing, the device shall be immediately sent for repair to an authorized service center.
- The use of excessive force during work shall be avoided.



### **WARNING:**

- During the work with the pipet controller general safety regulations regarding risks related with laboratory work should be observed. Protective clothing, goggles, and gloves should be worn.
- The pipet controller shall be used only for measuring liquids in conditions specified by the manufacturer, which are limited due to the chemical and mechanical resistance of the device, as well as the user safety.
- The information and instructions provided by the manufacturers of the reagents must be observed.
- Read the Instruction Manual before use. Improper handling of a Li-ion battery may result in the battery catching fire, overheating, or leaking.
- Do not puncture, crush, bend or modify the battery in any way.
- Never attempt to open or modify the battery. This can cause a fire or explosion.
- Do not expose the battery to excessive heat, direct sunlight, or fire.
- If there are signs of damage, swelling or leakage, stop use and contact the manufacturer immediately. Dispose the battery immediately according to rules in your country.
- Do not short circuit the (+) and (-) terminals with other metals.

**NOTE:** The pipet controller is equipped with liquid vapors exhaustion system which protects against corrosion to ensure long instrument life.

### 3.0 Limitations of Use


- The pipet controller shall not be used for measuring substances with vapors of which damage the following plastics: PP, SI, EPDM, PVDF, PTFE, PA.
- The pipet controller shall not be used in an environment where explosion risk is present.
- Flammable liquids shall not be measured – in particular, substances with flashpoint below 0°C (ether, acetone).
- The pipet controller shall not be used for drawing acids with a concentration above 1 mol/L.
- The pipet controller shall not be used for drawing solutions with a temperature above 50°C.
- The pipet controller may work in temperature range from +10°C to +35°C.

The pipet controller is suitable for general laboratory use only. It must be used only by personnel who know the health risks associated with the substances that are normally used with this instrument.

### 4.0 Switching On

The pipet controller is switched on by pressing any of the buttons. The display will show the selected device speed level, dispense mode and the battery level indicator.

Examples of the display indications:

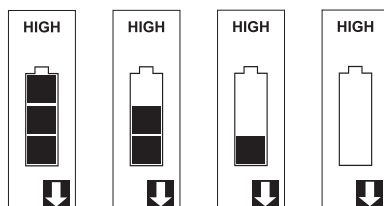

The battery is discharged and require charging if the indicator does not display “bars”. When the battery is fully charged, the indicator displays three “bars”.


- The pipet controller switches off automatically if not used for 5 minutes.
- The pipet controller may be charged only with the original charger.
- The mains voltage shall conform with the specification on the charger.
- Charging shall be done in accordance with Section 8 of this Instruction Manual.

## 5.0 Aspirating and Dispensing Liquids

### Setting device speed

The speed level (LOW or HIGH) should be set before aspirating is started. To change and preview the selected level/speed value press the Speed button.

Example: Selected level/speed value, HIGH 2 (H2).

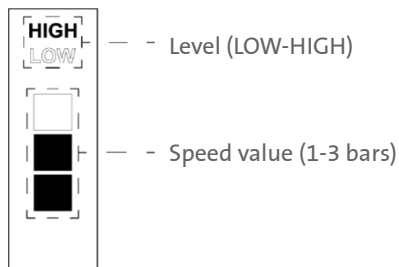

The liquids can be aspirated with two speed levels with 1 of 3 predefined values assigned for each of them:

- LOW: L1, L2, L3 (recommended for pipets of volumes up to 5 mL)
- HIGH: H1, H2, H3 (recommended for pipets of volumes greater than 5 mL)

**NOTE:** H3>H2>H1>L3>L2>L1. Step between the time of aspirating 25 mL of distilled water using the 25 mL Corning® Stripette® pipet on adjacent speed values is less than 1 s.

To change the speed value assigned to the current speed level:

- Press the Speed button for 3 seconds. The display will start flickering.
- Press the Speed button until the display shows the desired speed (1 to 3 bars).
- Press the Speed button again for 3 seconds. The display will flicker faster confirming new speed has been set.

If the Speed button is not pressed for 10 seconds or one of operating buttons is pressed, the speed reverts to the previous value.


## Attaching a pipet

**CAUTION:** Before attaching a pipet, check whether the pipet is not damaged, has no dents or sharp edges in the gripping part. Check whether the gripping part is dry.

The pipet shall be gripped as close to the upper end as possible and carefully inserted into the pipet holder until resistance is noticed.



**WARNING:** Do not apply excessive force so as not to damage thin pipets and avoid risk of injury.

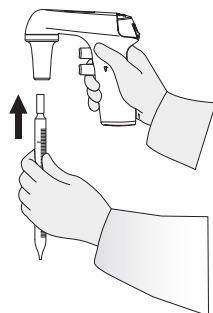

A pipet that has been correctly attached and sealed in the holder should not tilt to the sides.


After attaching a pipet, hold the pipet controller in a vertical position. It is not recommended to leave the device with a pipet attached for a longer period, for example overnight or over a weekend.

**CAUTION:** Do not put aside the pipet controller if there is liquid in the pipet.

## Filling the pipet

Holding the pipet controller in a vertical position, immerse the pipet end in the liquid to be drawn up, and press the aspiration button gently. The speed depends on how deep the aspirating button is pressed. The deeper the button is pressed the faster the liquid is aspirated into the pipet.

It is recommended to draw a slightly greater liquid volume than required (due to meniscus above the required volume mark), adjusting the aspiration speed, so as not to overfill the pipet.

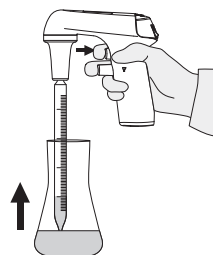


## Setting the volume

After the pipet is filled, dry the outside surface with absorbent paper that does not leave impurities. Then set the required liquid volume precisely. Pressing the dispense button gently, dispense the excessive liquid from the pipet until the meniscus of the liquid aligns exactly with the required volume mark on the pipet.

## Emptying the pipet

Holding the vessel in an inclined position, place the pipet end in contact with the vessel wall and press the dispense button gently. The dispensing intensity may be adjusted depending on how deep the dispense button has been pressed. The deeper the button is pressed the faster the outflow of liquid from the pipet.

The pipet controller has two dispense modes. The dispense mode is selected by successive pressing the Mode button until the display shows the correct mode.

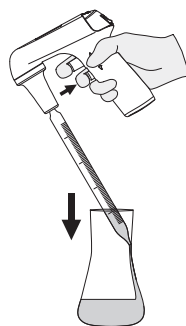


• **Gravity mode** 

Dispensing is effected in gravity mode, which means that the liquid flows out of the pipet by its own weight.

• **Blow out mode** 

Dispensing is effected in gravity mode, however, when the dispense button is pressed to the middle position, the pump is started and fast emptying of the pipet with a blow out is effected.

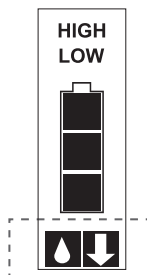

**CAUTION:** During gravimetric dispensing the pipet is not completely emptied due to the characteristics of pipets used with the pipet controller.

## 6.0 Troubleshooting

If during your work the pipet controller is not functioning correctly, check the cause and correct the fault.

Problem	Possible Cause	Action
The pipet falls out (the holding force of the pipet is too small), or tilts to the side too much.	The pipet holder is dirty or wet.	Take out the pipet holder, and clean, wash, and dry it.
	The pipet holder is damaged.	Replace the pipet holder with a new one.
The pump is working, but the pipet controller aid does not draw liquid or draws liquid very slowly.	The filter is dirty.	Take out the pipet holder, take out the filter; if it is dirty, replace it with a new one.
	The pipet holder and/or the connector gasket is damaged.	Replace the mechanically damaged elements with new ones.
	The pipet is damaged.	Check the pipet for damage (cracks, dents); if present, replace the pipet with a new one.
Liquid leaks from the pipet (the aspiration and the dispense buttons are not pressed).	The pipet is inserted incorrectly.	Check whether the pipet has been correctly inserted in the pipet holder.
	The pipet holder, the filter, or the connector gasket is installed incorrectly.	Check whether all parts are present and correctly installed.
	The pipet holder and/or the connector gasket is damaged.	Replace the mechanically damaged elements with new ones.


If the above actions do not help and service is required, contact Corning Customer Service. Before sending for service, the pipet controller should be cleaned and decontaminated. Written details including the precise specification of solutions used and the type of laboratory in which the device was used, should be sent with the product.

## 7.0 Replacing the Filter

**CAUTION:** The work safety instructions given in Section 2 shall be observed when disassembling the pipet controller.

The filter replacement is necessary, if drawing efficiency deterioration is observed. The direct reason may be a dirty filter after a long period of use.

In order to replace the filter:

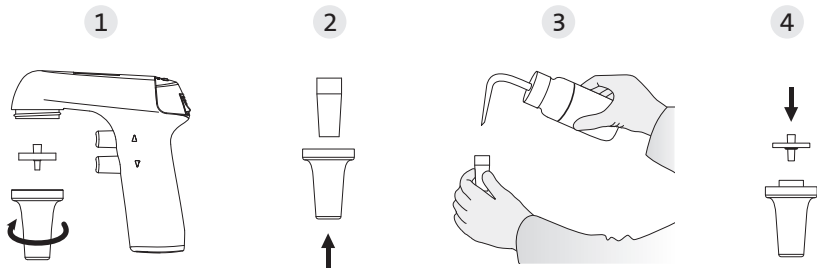


1. Remove the pipet. Unscrew the nose piece and remove the membrane filter.
2. Remove the pipet holder.
3. Rinse the holder using a wash bottle. Blow liquid out of the holder and set it aside until it is completely dry.
4. Install a new membrane filter and assemble the device in reverse order.

## 8.0 Charging the Battery

The Corning® Stripettor® Pro Pipet Controller (Cat. Nos. 4999, 4900) comes with a universal charger and a set of adapters in different versions: EU, USA, GBR, KOR, and AUS. Choose your country's adapter and connect to the housing.

EU and KOR adapters look very similar and are distinguished with dedicated stickers.

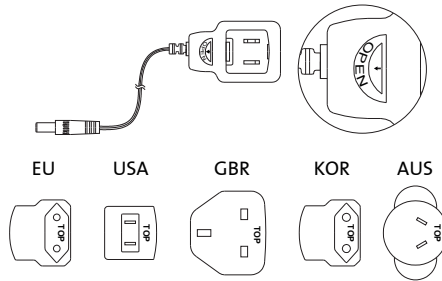

To mount the adapter, it should be inserted into the slots of the housing in the direction of the arrow, until you hear a click.

To remove or change the adapter, simply press the Open button in the direction of the arrow, hold the button down, remove the adapter in the direction of the arrow.

**CAUTION:** The pipet controller may be charged only with the original charger. The mains voltage shall conform with the specification on the charger (Input: AC 100-240V, 50/60 Hz). Using chargers other than the original one may damage the battery.

The pipet controller is delivered with 1 Li-ion battery.

### Charging



#### WARNING:

- Stop charging the battery if charging is not completed within the specified time.
- Stop using the battery if abnormal heat, odor, discoloration, deformation or abnormal condition is detected during use, charge, or storage. Dispose the battery immediately according to rules in your country.

#### WARNING: In order to prolong the life span of the rechargeable battery, the following rules should be followed:

- Before the pipet controller is activated for the first time, the battery should be charged.
- If the pipet controller starts to indicate low battery level during work, connect it to the power supply to continue working.
- Do not leave the pipet controller fully discharged for a long period of time (more than 1 week).
- In case of a planned break longer than 1 month in using the device, it is recommended to charge the battery fully.

1. Charging temperature: 10°C to 35°C.
2. Charging the battery is carried out through a charger (power supply) by direct connection to the main power. Battery charging is indicated by successive lighting of “bars”.
3. Full charging time: approx. 3 hours.
4. The battery is charged when all 3 “bars” are displayed simultaneously.


When the battery is fully charged, the charging circuit disconnects automatically. If used correctly, the service life of the battery should be at least 500 full charging cycles before it's capacity drops below 80% of rated capacity. If the battery shows signs of wear and not holding the charge long enough, replace the battery. If all instructions of the manufacturer are followed, it is not possible to overcharge the battery.

### Charging on the charging stand

1. Put the pipet controller on the stand.
2. Plug the charger into the stand outlet.
3. Connect the charger to the adapter and to the main power.

## 9.0 Replacing the Battery

If necessary, the battery can be replaced easily.

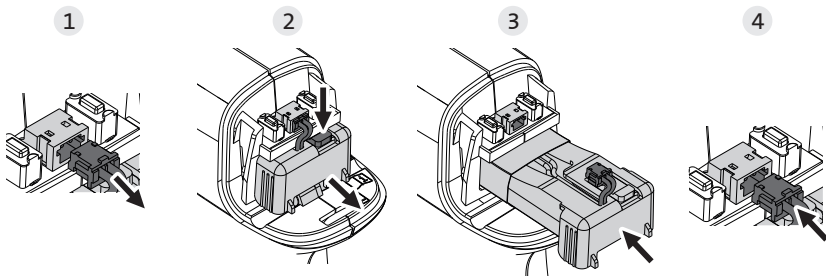

1. Take off the battery cover. Unplug the battery cable from the socket.
2. Remove the battery set from the battery compartment.
3. Slide in the new battery set.
4. Plug in the battery cable to the socket.
5. Close the battery cover.

## 10.0 Maintenance

### Cleaning

The pipet controller does not require any maintenance. Its external parts may be cleaned with a swab moistened with isopropyl alcohol.

The nose piece and the pipet holder may be autoclaved at 121°C for 20 minutes.

After autoclaving, dry the pipet holder. The filter included in the set may be sterilized by autoclaving at 121°C for not more than 15 minutes.

### Ultraviolet (UV) sterilization

The outer body of the pipet controller is UV-resistant, which was confirmed by many tests. The recommended distance from the radiation source to exposed element should be not less than 50 cm.


Prolonged and very intense UV exposure can cause discoloration of pipet controller parts, without affecting its performance.

### Storage

The pipet controller should be stored in a dry place. The allowable storage temperature is -20°C to +45°C.


During breaks in the work the pipet controller can be stored on the wall hanger or charging stand. The Corning® Stripettor® Pro charging stand is a device intended to conveniently charge and store a pipet controller when not in use. Three positions to store the pipet controller, depending on the user needs, are also visible as icons printed on the stand.

**CAUTION:** Do not store the pipet controller with a filled pipet.

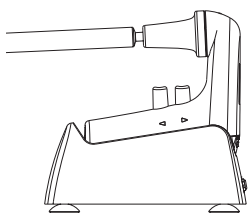


#### Break position

Convenient while working, helps to avoid contamination.

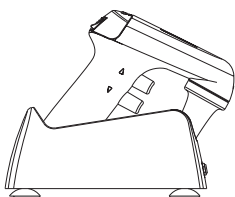

#### Rest position

Ideal way to store a pipet controller after use.

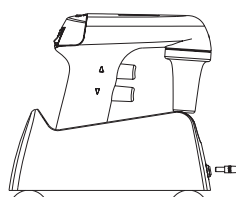

#### Charging position

Used while pipet controller is in charging process.

## 11.0 Components

The pipet controller set is supplied with the following components:

- Li-ion battery set
- PTFE filter 0.2 µm
- Wall hanger
- Charging stand, 3-positions (included with Cat. No. 4999)
- Universal charger with set of adapters
- Quick Start Guide (CLS-AN-818DOC)

## 12.0 Spare Parts

Cat. No.	Description	Qty/Pk
4997	Set of 4 colored nose pieces (orange, yellow, green, red)	4
4993	Silicone pipet holder	1
4991	PTFE filter 0.2 $\mu\text{m}$	5
4992	PTFE filter 0.45 $\mu\text{m}$	5
4947	Li-ion battery set	1
4948	Battery cover	1
4998	Charging stand, 3-positions	1
4996	Universal power supply, 9V with set of adapters (EU, USA, GBR, KOR, and AUS)	1

## 13.0 Limited Warranty

Corning Incorporated (Corning) warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. CORNING DISCLAIMS ALL OTHER WARRANTIES WHETHER EXPRESSED OR IMPLIED, INCLUDING ANY IMPLIED WARRANTIES OF MERCHANTABILITY OR OF FITNESS FOR A PARTICULAR PURPOSE. Corning's sole obligation shall be to repair or replace, at its option, any product or part thereof that proves defective in material or workmanship within the warranty period, provided the purchaser notifies Corning of any such defect. Corning is not liable for any incidental or consequential damages, commercial loss or any other damages from the use of this product.

This warranty is valid only if the product is used for its intended purpose and within the guidelines specified in the supplied instruction manual. This warranty does not cover damage caused by accident, neglect, misuse, improper service, natural forces or other causes not arising from defects in original material or workmanship. This warranty does not cover batteries or damage to paint or finish. Claims for transit damage should be filed with the transportation carrier.

In the event this product fails within the specified period of time because of a defect in material or workmanship, contact Corning's Customer Service at the following numbers: USA: 1.800.492.1110; Canada: 1.978.442.2200. For other regions of the world, visit [www.corning.com/lifesciences](http://www.corning.com/lifesciences) or see the included instruction manual for a list of Worldwide Support Offices.

Corning's Customer Service team will help arrange local service where available or coordinate a return authorization number and shipping instructions. Products received without proper authorization will be returned. All items returned for service should be sent postage prepaid in the original packaging or other suitable carton, padded to avoid damage. Corning will not be responsible for damage incurred by improper packaging. Corning may elect for onsite service for larger equipment. Some states do not allow limitation on the length of implied warranties or the exclusion or limitation of incidental or consequential damages. This warranty gives you specific legal rights. You may have other rights which vary from state to state.

No individual may accept for, or on behalf of Corning, any other obligation of liability, or extend the period of this warranty.

For your reference, make a note of the serial and model number, date of purchase, and supplier here.

Serial No. \_\_\_\_\_ Date Purchased \_\_\_\_\_

Model No. \_\_\_\_\_

Supplier \_\_\_\_\_

## 14.0 Equipment Disposal



According to Directive 2012/19/EU of the European Parliament and of the Council of 4 July 2012 on waste electrical and electronic equipment (WEEE), this product is marked with the crossed-out wheeled bin and must not be disposed of with domestic waste.

Consequently, the buyer shall follow the instructions for reuse and recycling of waste electronic and electrical equipment (WEEE) provided with the products and available at [www.corning.com/weee](http://www.corning.com/weee).

**How to Purchase:** For specific availability in your region and purchasing options, terms and conditions of sale, customer/product support, and certificates, visit [www.corning.com/how-to-buy](http://www.corning.com/how-to-buy).

**Warranty/Disclaimer:** Unless otherwise specified, all products are for research use or general laboratory use only.\* Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. These products are not intended to mitigate the presence of microorganisms on surfaces or in the environment, where such organisms can be deleterious to humans or the environment. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications. \*For a listing of US medical devices, regulatory classifications or specific information on claims, visit [www.corning.com/resources](http://www.corning.com/resources).

*Corning's products are not specifically designed and tested for diagnostic testing. Many Corning products, though not specific for diagnostic testing, can be used in the workflow and preparation of the test at the customers discretion. Customers may use these products to support their claims. We cannot make any claims or statements that our products are approved for diagnostic testing either directly or indirectly. The customer is responsible for any testing, validation, and/or regulatory submissions that may be required to support the safety and efficacy of their intended application.*

# CORNING

**Corning Incorporated**  
*Life Sciences*

[www.corning.com/lifesciences](http://www.corning.com/lifesciences)

## **NORTH AMERICA**

t 800.492.1110

t 978.442.2200

## **ASIA/PACIFIC**

### **Australia/New Zealand**

t 61 427286832

### **Chinese Mainland**

t 86 21 3338 4338

### **India**

t 91 124 4604000

### **Japan**

t 81 3-3586 1996

### **Korea**

t 82 2-796-9500

### **Singapore**

t 65 6572-9740

### **Taiwan**

t 886 2-2716-0338

## **EUROPE**

[CSEurope@corning.com](mailto:CSEurope@corning.com)

### **France**

t 0800 916 882

### **Germany**

t 0800 101 1153

### **The Netherlands**

t 020 655 79 28

### **United Kingdom**

t 0800 376 8660

### **All Other European Countries**

t +31 (0) 206 59 60 51

## **LATIN AMERICA**

[grupoLA@corning.com](mailto:grupoLA@corning.com)

### **Brazil**

t 55 (11) 3089-7400

### **Mexico**

t (52-81) 8158-8400