


INSTRUCTIONS

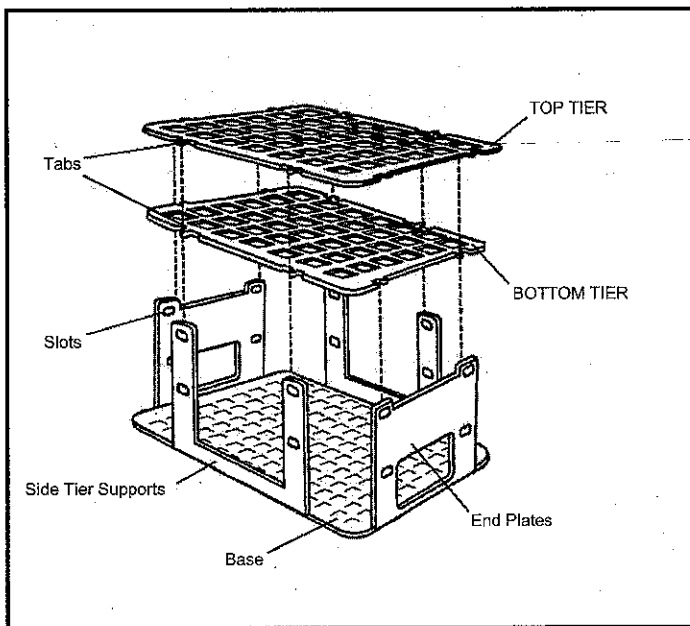

- CAN BE ERECTED IN WATER
- AUTOCLAVABLE
- STACKABLE
- EASY TO ASSEMBLE
AND TAKE APART
- TUBES EASY TO IDENTIFY
- PLATES PROVIDED
FOR LABELING

1. Place the base on a flat work surface with the word "BOTTOM" facing down.
2. Bend up one end plate. Snap the bottom tier end tabs into the lower set of slots in the end plate. Be sure the bottom tier imprinted numbers face up.
3. Position the top tier so that imprinted numbers face up and align with the numbers on the bottom tier. Snap the top tier end tabs into the upper set of slots in the end plate.
4. Bend up the opposing end plate and snap the bottom tier and then top tier end tabs into their matching slots in the end plate.
5. Bend up one side tier support and snap the bottom tier and then top tier side tabs into their matching slots in the side tier supports.
6. Bend up the remaining side tier support and snap the bottom tier and then top tier side tabs into their matching slots in the side tier supports.