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Operation Manual

- Switch to clock mode or timer mode function : Press [タイマー/時計] button
- Clock setting
 1. Press [タイマー/時計] button to enter clock mode.
 2. Keep pressing [タイマー/時計] for 3 seconds until sounds [PiPi], then the digit in hour position will blink.
 3. Press the [10 秒] button to select AM, PM, 24-hour display.
 4. Press [10 分] button and [1 分] button to set hour value.
 5. Press [タイマー/時計] button to enter the minute setting, then the digit in minute position will blink.
 6. Press [10 分] button and [1 分] button to set minute value.
 7. Press [タイマー/時計] button again to finish the clock setting.
- Counting down
 1. Press [タイマー/時計] button to enter timer mode.
 2. Press [リセット] button to display [00MOOS].
 3. Press [10 分] [1 分] [10 秒] button to set the timer.
Example) setting counting down time at 21 min30sec : press [10 分] button 2 times → press [1 分] button 1 time → press [10 秒] button 3 times.
 4. Press [スタート/ストップ] to start counting down. After time is up, the timer sounds for 15 seconds.
 5. Press any key to stop the sound.
(Repeat function) : Returning to previous count down time setting after stopping alarm.
 6. Press [スタート/ストップ] button to pause counting down during counting down. And press [スタート/ストップ] button again to restart counting down from the paused time.
 7. Press [タイマー/時計] button to switch to clock mode during counting down. After counting down to zero, this timer sounds in clock mode as well.
- Counting up
 1. Press [タイマー/時計] button to enter timer mode.
 2. Press [リセット] button to display [00MOOS].
 3. Press [スタート/ストップ] button to start counting up.
 4. Press [スタート/ストップ] button to stop counting up. Press [スタート/ストップ] button again to restart counting up from the paused time.
 5. Press [リセット] button to reset the counting and show 00MOOS.
※After the counting up to [99M59S], the timer will return to [00MOOS] display and re-counting up from zero.

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