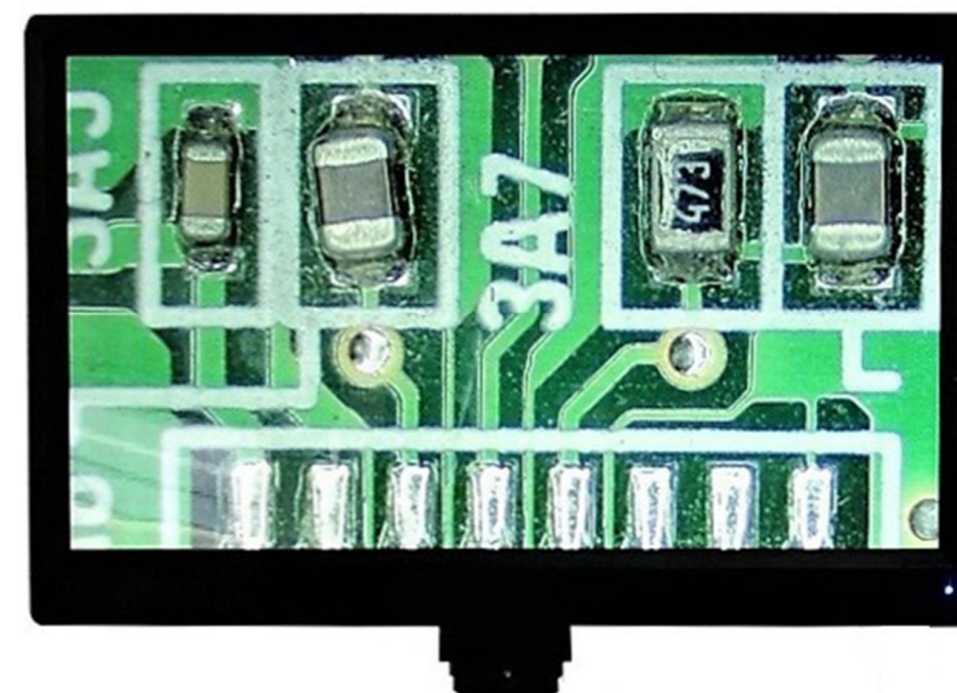
INSTRUCTION MANUAL

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XR9027 / DP300
LCD Screen Camera
LCDカメラ

Specifications (仕様)

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T : Text annotation: Click the button to bring up the checkbox for character annotation.

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 : Feature point: After clicking the button, click any point in the interface to determine the starting point, and click a second time to confirm the direction and position of the arrow.

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 : Color: Click this button to select the color of the line, eight colors available.

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 : Line width: Click this button to select the thickness of the line, a total of 4 sizes optional.

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 : Clear screen: When the mouse clicks on this button, all items on the screen are deleted.

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 : Undo: Mouse click on this icon to undo the previous operation.

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 : Export: Click this button to save the file to the storage space.

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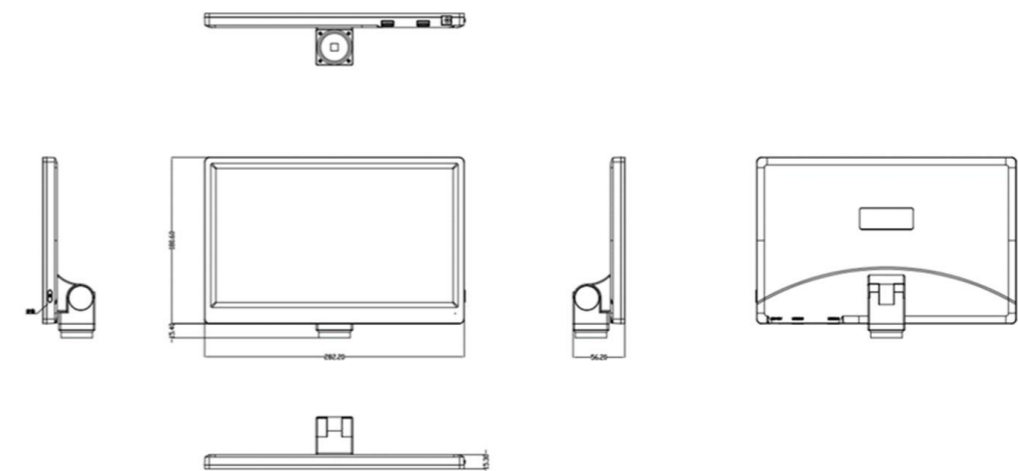
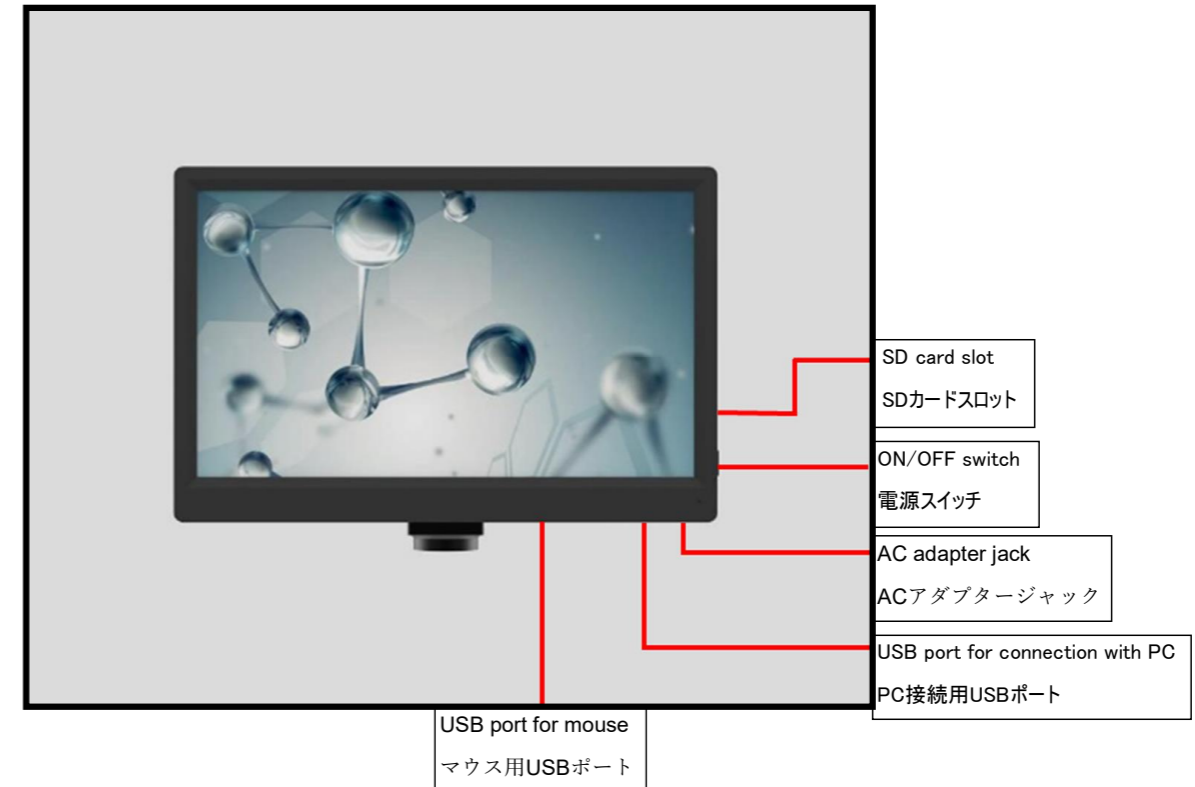
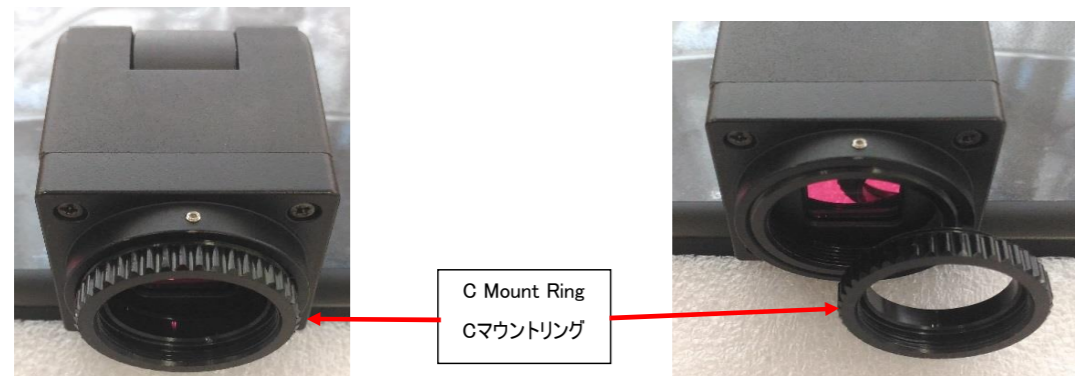
DP300 is provided with a C mount ring to be mounted on DSZV-44C, SPZT-50, SPZV-50 and SPZHT-135 microscopes (C mount). When mounting it on DSZV-44 microscopes (CS mount), unscrew the C mount from the camera.

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Take out an USB receiver from the mouse and plug the receiver into the USB port where a mouse is illustrated on the back of the device. No battery for the mouse is included.

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3. Power on the device (電源入力)

Connect an AC adapter with the device, and plug in the AC adapter. A power indicator lamp turns red, and press and hold an on/off switch until the power indicator lamp turns blue.

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When connecting the USB port to PC with a Type A - Type A USB cable, you can use it as a normal UVC camera.

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 : Radius fixed circle: after the mouse clicking the button, click any point on the interface to determine the center of the circle, the second click to determine the distance from the center of the circle. The measurement is displayed as the radius, area and circumference of the circle.

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 : Two-point fixed circle: after the mouse clicking the button, click any point on the interface to determine the point of the circle circumference. Click again to determine the circle, and the measurement is displayed as the circumference and area of the circle.

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 : Three points fixed circle: After clicking the button with the mouse, click any three points on the interface to determine the circumference and area of the circle.

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 : Concentric circle: After clicking the button with the mouse, click any point on the interface to determine the center of the circle, the second click to determine the first circle, and the third click to determine the second circle. The measurement is displayed as the distance between the two circles.

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 : Point-circle distance: after clicking the button, click any point in the interface to determine the center of the circle, the second click to determine the circle, the third click to determine the distance between the line and the center of the circle. The measurement is displayed as the distance from the point to the center of the circle.

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 : Circle centers distance: After clicking the button, click any point in the interface to determine the center of the circle, the second click to determine the first circle, the third click to confirm the center of the second circle, and the fourth click to confirm the second circle. The measurement is displayed as the distance between the centers of the two circles.

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 : Vertical line distance: after the mouse clicking the button, click any point on the interface to determine the starting point, the second click to determine the straight line, the third click to determine the vertical distance from the line, measuring the length of the vertical line (after confirming the vertical line segment, drag the length shown by the cursor to the appropriate location).

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 : Rectangle: After clicking the button with the mouse, click any point on the interface to determine the right angle position, click second time to determine the perimeter of the rectangle, measure the perimeter and area of the rectangle (after confirming the perimeter, drag the perimeter and area shown by the cursor to the appropriate location).

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 : Polygon: after clicking the button with the mouse, click any point on the interface to determine a corner. Each click increases a corner, and the last click should overlap with the first clicked point, measure the perimeter and area occupied by the polygon (after confirming OK, drag the perimeter and area shown by the cursor to the appropriate location).

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 : Angle measurement: after the mouse clicking the button, click any point on the interface to determine the starting point, the second click to determine the straight line, the third click to determine the straight line between the second point and measure the angle between the two straight lines (to confirm the angle, drag the angle displayed by the cursor to the appropriate location).

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 : Angle measurement 2): Click 2 points on one straight line and also click 2 points on another straight line to measure the angle of the two lines. The measured value is stuck on the place where you click after the measurement.

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 : Arc: after the mouse clicking the button, click any point on the interface to determine the starting point, the second click to determine the straight line, the third point to determine the straight line with the second point. measure the length of the arc by three points of arc.

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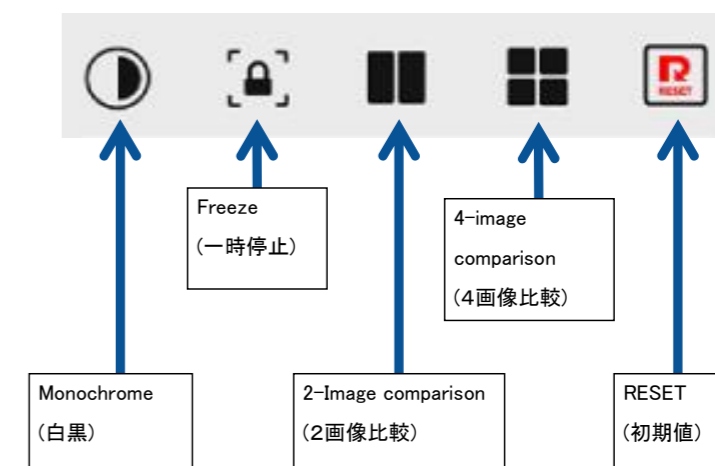
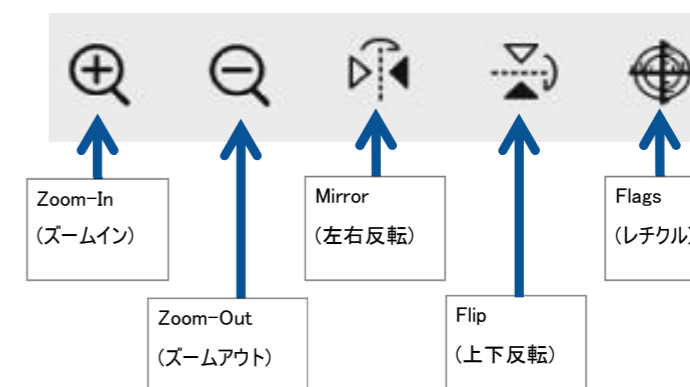
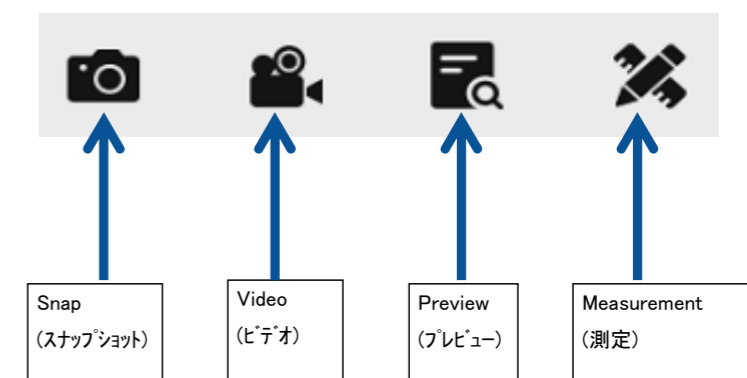
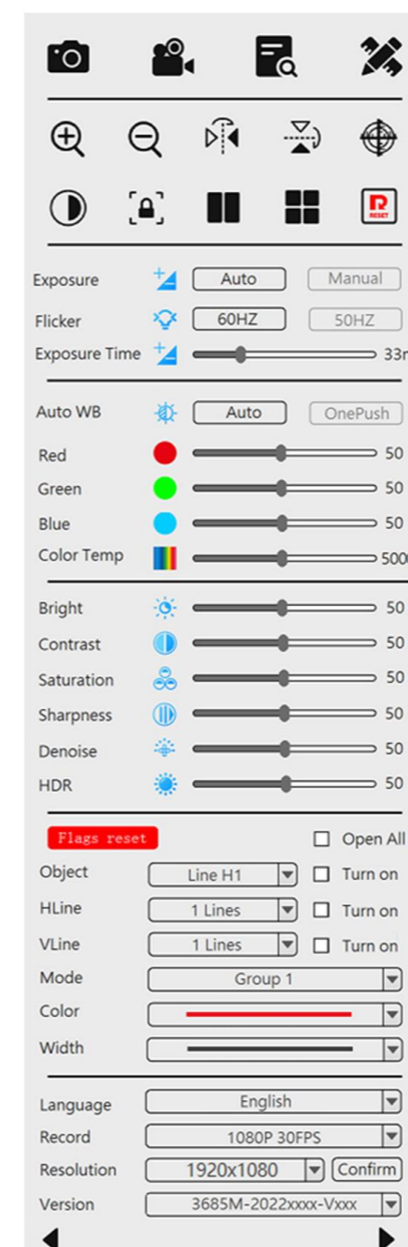
Connect a mouse with the device. When moving a cursor to the left (or right) side on the screen, the main menu will pop up as follow.

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Two-image comparison (2画面比較)

The saved image(s) and the live image are displayed when clicking this icon. Click a saved image to compare it with the live image.

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Four-image comparison (4画面比較)

Four live images are displayed when clicking this icon. One of the four live images is frozen by double-clicking and back to a live image by double-clicking again.

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Flicker (フリッカー)

Select the frequency in your area, 60HZ or 50HZ

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Exposure Time (露出時間)

Adjust the exposure time by dragging the slider.

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
Auto is to set all of the three colors, R,G,B at 50%. OnePush is to adjust white balance automatically on the basis of the white object displayed. When adjusting manually, drag the correspondent sliders.

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Measurement Tools (計測ツール)

Click the icon of the measurement function that you want to use, and after measurement, right-click to end the measurement.

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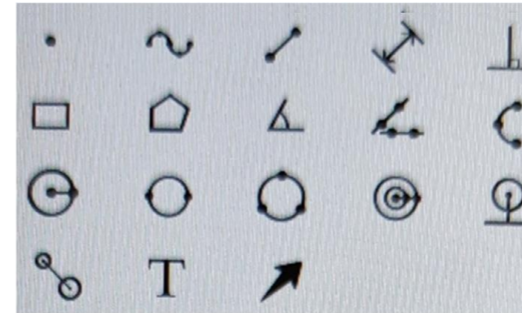 : Point Coordinate: After clicking the button with the mouse, click OK at any point in the interface to draw a point and display the coordinates of the point.

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 : Curve: after the mouse clicking the button, click any point on the interface to determine the starting point. The mouse passes through the path for the length, and the second click determines the end point and measures the length of the drawn line.

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 : Point line distance: After mouse clicking the button, click any point in the interface to determine the starting point, and click a second time to determine the distance between the starting point and the end point. Measure the length of the line (after drawing the line, then click the mouse at the appropriate location, the length of the line will be displayed).

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 : Lines distance: after clicking the button, click any point in the interface to determine the starting point, the second click to determine the straight line, the third click to determine the position of the parallel straight line. Measure the distance between the two parallel lines (after drawing the line, drag the length that the cursor shows to the appropriate position).

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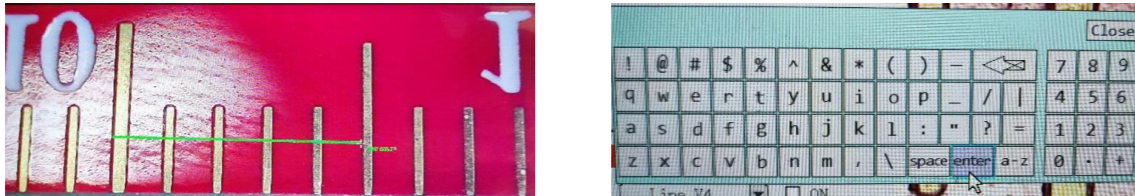
After displaying a calibration scale and focusing on it, click the line calibration icon to make a calibration. Click two points to draw a calibration line on the basis of the increments of the calibration scale displayed. Input the calibration name and the calibration length with the key board which is popped up when clicking each column, and also select the unit and the accuracy from drop down lists. Make sure to save the calibration by clicking the save icon.

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When deleting the calibration set, click the delete icon.

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After calibration, make sure to save it by clicking.

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Image Adjustment (画質調整)

These sliders allow you to adjust image parameters.

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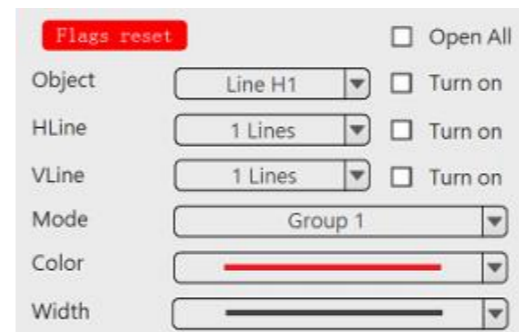
You can set 8 kinds of grid lines. The grid lines can be moved by dragging when necessary.

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To select the line which you want to change its color and/or width.

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To select quantity of horizontal line.

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Mode(モード)

To set and save 8 kinds of grid lines in Group 1 to 8.

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Language

Record

You can select a language from English, Chinese, Taiwanese, Japanese and Spanish. There are two resolutions selectable, 1080P or 720P.

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Version

The main menu moves to the right or left side where the triangle icon is clicked.

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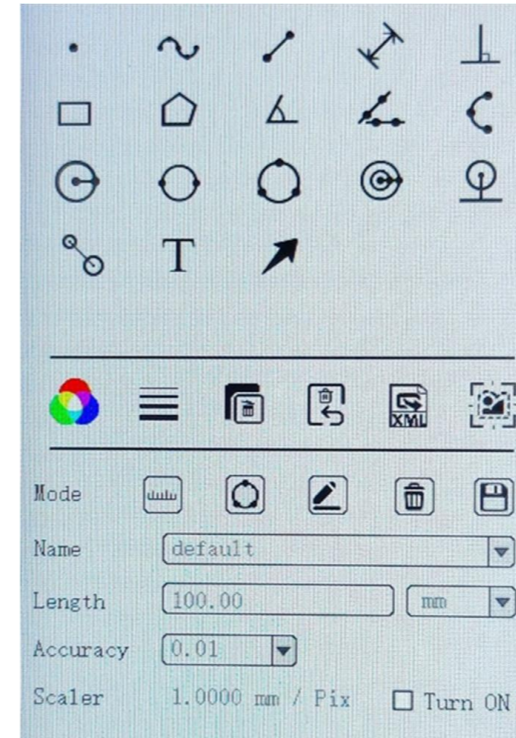
Before measurement, you need to make calibration according to the magnification of the microscope that you use.

*) Note: Measurement function can not be used without calibration.

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